




7585

Material **Nitrile** LENGTH **12 in. / 300mm**

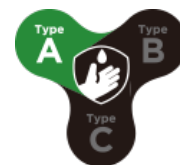

CHEMICAL PERMEATION

CHEMICAL NAME	CAS NUMBER	BDT
		TTL EN374
Formaldehyde 37%	50-00-0	>120
2-Hydroxypropionic acid 85%	50-21-5	>480
Carbon Tet	56-23-5	6-10
1,2-Propanediol	57-55-6	>480
Diethyl Ether	60-29-7	1-5
Aminobenzene	62-53-3	6-10
Ethanol	64-17-5	>30
Formic Acid 90%	64-18-6	1-5
Acetic Acid 99%	64-19-7	6-10
Acetic Acid 50%	64-19-7	>60
Acetic Acid 84%	64-19-7	>10
2-Propanol	67-63-0	>120
2-Propanone	67-64-1	<1
Chloroform	67-66-3	<1
Dimethylsulfoxide (DMSO)	67-68-5	>10
Dimethyl Formamide	68-12-2	1-5
n-Propanol	71-23-8	>60
Butanol	71-36-3	>60
Alcohol, Amyl	71-41-0	>60
Benzene	71-43-2	1-5
Iodide, Methyl	74-88-4	<1
ETHYLAMINE 70%	75-04-7	<1
Acetonitrile	75-05-8	1-5
Acetaldehyde	75-07-0	<1
Chloride, Methylene	75-09-2	<1
Carbon Disulfide	75-15-0	<1

Bromoform	75-25-2	<1
1,1-Dichloroethene	75-35-4	1-5
Acetyl Chloride	75-36-5	1-5
Nitromethane	75-52-5	<1
1,2-Epoxypropane	75-56-9	<1
Tetramethylammonium Hydroxide 25%	75-59-2	>480
Dimethyl Sulfate	77-78-1	>10
Citric Acid 99%	77-92-9	>480
Citric Acid 50%	77-92-9	>480
Citric Acid 75%	77-92-9	>480
Citric Acid 30%	77-92-9	>480
Dichloropropane, 1,2-	78-87-5	1-5
2-Butanone	78-93-3	1-5
Ethylene, Trichloride	79-01-6	1-5
2-Propeneamide 98%	79-06-1	>480
Acetate, Methyl	79-20-9	<1
PERACETIC ACID 39%	79-21-0	1-5
Nitro Propane	79-46-9	<1
Methacrylate, Methyl	80-62-6	1-5
DBP	84-74-2	>120
Vinyl Pyrrolidinone	88-12-0	1-5
Dichlorobenzene O-	95-50-1	1-5
2-Aminotoluene	95-53-4	1-5
Pseudocumene	95-63-6	1-5
Butanone Oxime	96-29-7	>60
2-Ethylbutyl alcohol	97-95-0	1-5
2-Furaldehyde	98-01-1	1-5
Butyl Toluene, p-Tert	98-51-1	>30
(1-Methylethyl)benzene	98-82-8	1-5
Nitrobenzene	98-95-3	1-5
Benzene, Ethyl	100-41-4	1-5
Benzene, Vinyl	100-42-5	1-5
Alcohol, Benzyl	100-51-6	1-5
Benzaldehyde	100-52-7	<1
2,2',2''-Nitrilotriethanol	102-71-6	>30
1,2-Dichloroethane	107-06-2	1-5
Acrylonitrile	107-13-1	1-5

1,2-Diaminoethane 99%	107-15-3	1-5
2-Propen-1-ol	107-18-6	6-10
1,2-Ethanediol	107-21-1	>480
Methyl Propyl Ketone	107-87-9	1-5
1-methoxy-2-propanol	107-98-2	>10
Acetate, Vinyl	108-05-4	<1
2-Pentanone, Methyl-	108-10-1	1-5
Acetate, Isopropyl	108-21-4	1-5
PROPYLENE GLYCOL MONOMETHYL ETHER ACETATE	108-65-6	6-10
2,6-Dimethyl-4-Heptanone	108-83-8	>10
Benzene, Methyl	108-88-3	1-5
Benzene Chloride	108-90-7	1-5
Cyclohexanol	108-93-0	>120
Cyclohexanone	108-94-1	1-5
Carbolic Acid(s) 100%	108-95-2	>30
Carbolic Acid 10%	108-95-2	6-10
Carbolic Acid 89%	108-95-2	6-10
Dimethyl Propaneamide, N,N'-	109-55-7	1-5
Acetate, Propyl	109-60-4	1-5
Pentane	109-66-0	>120
1-Aminobutane	109-73-9	<1
Diethylamine	109-89-7	1-5
Diethylene Oxide	109-99-9	1-5
Hexane	110-54-3	>240
Cyclohexane	110-82-7	>240
PYRIDINE	110-86-1	<1
Diethylene oximide	110-91-8	<1
1,5-Pentanedial 50%	111-30-8	>60
2,2-iminodiethanol	111-42-2	>30
n-Octane	111-65-9	>480
2-Butoxyethanol	111-76-2	>30
n-Octanol	111-87-5	>30
Oleic Acid 98%	112-80-1	>480
1,2,4-Trichlorobenzene	120-82-1	1-5
TRIETHYLAMINE	121-44-8	>30
4-Hydroxy-4-methyl-2-pentanone	123-42-2	1-5
3-Methyl-1-butanol	123-51-3	>60

Butyl Acetate	123-86-4	1-5
1,4-Dioxane	123-91-1	1-5
3-Methylbutyl Ethanoate	123-92-2	<1
PERC	127-18-4	1-5
Dimethylacetamide N,N-	127-19-5	1-5
Butyl Acrylate	141-32-2	1-5
2-Aminoethanol	141-43-5	1-5
Ethyl Acetate	141-78-6	1-5
Heptane	142-82-5	>480
Butoxytriglycol	143-22-6	>10
OXALIC ACID (s) 99%	144-62-7	>480
Calcium Carbonate (s) 99%	471-34-1	>480
Trimethyl Phosphate	512-56-1	6-10
2,2,4-Trimethyl Pentane	540-84-1	>240
3-Methyl-2-Butanone	563-80-4	1-5
Butyl Ethylene	592-41-6	>10
Amyl Acetate	628-63-7	1-5
Methyl Pyrrolidone, N-	872-50-4	6-10
2-Bromoethyl Acetate	927-68-4	1-5
Hexamethyldisilazane	999-97-3	>480
3,8-Diamino-5-ethyl-6-phenylphenanthridinium bromide 1%	1239-45-8	>480
3,8-Diamino-5-ethyl-6-phenylphenanthridinium bromide 5%	1239-45-8	>480
3,8-Diamino-5-ethyl-6-phenylphenanthridinium bromide 95%	1239-45-8	>480
Calcium Hydroxide (s) 95%	1305-62-0	>480
Caustic Potash 45%	1310-58-3	>480
Caustic Potash 10%	1310-58-3	>480
Caustic Potash 99%	1310-58-3	>480
Caustic Soda 50%	1310-73-2	>480
Caustic Soda 10%	1310-73-2	>480
Caustic Soda 40%	1310-73-2	>480
Caustic Soda 98%	1310-73-2	>480
Cresols	1319-77-3	1-5
Divinyl Benzene	1321-74-0	1-5
dimethyl benzene	1330-20-7	1-5
Chromic Acid Solution 50%	1333-82-0	<1
Ammonia Solution 29%	1336-21-6	>30
Ammonia Solution 25%	1336-21-6	>60

Gallotannin 95%	1401-55-4	>480
Methyl-Tert-Butyl Ether	1634-04-4	>10
Butoxypropanol	5131-66-8	>60
D-Limonene	5989-27-5	6-10
Hydrochloric Acid 37%	7647-01-0	>240
Muriatic Acid 20%	7647-01-0	>480
Muriatic Acid 32%	7647-01-0	>480
Hydrochloric Acid 10%	7647-01-0	>480
Phosphoric Acid 85%	7664-38-2	>480
Phosphoric Acid 50%	7664-38-2	>480
Phosphoric Acid 10%	7664-38-2	>480
Hydrofluoric Acid 30%	7664-39-3	>60
Hydrofluoric Acid 40%	7664-39-3	>10
Hydrofluoric Acid 48%	7664-39-3	>10
Battery Acid 47%	7664-93-9	>480
Battery Acid 10%	7664-93-9	>480
Battery Acid 25%	7664-93-9	>480
Battery Acid 96%	7664-93-9	>30
Battery Acid 50%	7664-93-9	>480
Bleach: Sodium Hypochlorite 6%	7681-52-9	>480
Bleach: Sodium Hypochlorite 12%	7681-52-9	>480
Nitric Acid 50%	7697-37-2	>30
Nitric Acid 10%	7697-37-2	>480
Nitric Acid 65%	7697-37-2	1-5
Nitric Acid 70%	7697-37-2	1-5
Hydrogen Peroxide 30%	7722-84-1	>480
Hydroxylamine 50%	7803-49-8	>480
Fir Oil	8006-64-2	>240
Oleum (20% Free SO3)	8014-95-7	<1
Ligroin	8032-32-4	>10
Dry Cleaning Mineral Spirits	8052-41-3	>480
Hydrobromic Acid 48%	10035-10-6	>480
Tetrachloropropene	10436-39-2	1-5
Ammonium Fluroide 40%	12125-01-8	>480
2-Chloro-2-Oxoethyl Acetate	13831-31-7	1-5
Pentachloropropane	23153-23-3	6-10
Antimony Tributyratate 95%	53856-17-0	>480

Dry cleaning safety solvent	64475-85-0	>480
Kerosene (hydrodesulfurized)	64742-81-0	>480
Diesel Oil	68334-30-5	>480
Diesel Fuel #2	68476-34-6	>480

BDT=BREAKTHROUGH DETECTION TIME

THE LEVEL (0 TO 6) INDICATES THE TIME REQUIRED FOR DIFFERENT CHEMICALS TO PERMEATE THROUGH THE GLOVE.

TTL : TOTAL IMMERSION CHEMICAL PERMEATION BREAKTHROUGH TIME.

INT : INTERMITTENT CONTACT CHEMICAL PERMEATION BREAKTHROUGH TIME, ONE MINUTE IMMERSION OUT OF EVERY TEN, REPEATEDLY.

Warranty Limitations and Disclaimer Use

This information is provided solely as a convenience to help you evaluate our gloves in the end-user's particular application. It is the responsibility of the purchaser and/or user to determine the level of toxicity of the materials to be handled and to select the proper glove suitable for a particular application. The information provided reflects laboratory performance of gloves under carefully controlled conditions. SHOWA makes no guarantee of results and assumes no obligation or liability in connection with this information.